FermentoFlash

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The device for the automatic analysis of beer

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Measurement principle:

The beer sample (10 ml) is sucked into the measuring cells by means of a pump. Alcoholic content, extract and density are measured by using thermal measuring effects. Derived constituents as original wort, apparent extract and osmotic pressure also will be determined.

Calibration:

The device may be calibrated by using a corresponding reference beer. You can store 18 different sets of calibration data (for example Pils, Strong beer, Non alcoholic beer). Then you can change from one product to another without having to undertake a new calibration.

Interfaces:

Parallel interface / recording printer: The device has a parallel interface for connecting up a normal commercial printer. A recording printer is included in the scope of delivery.

Serial interface

A PC can be connected to the serial interface. The software that is necessary for this is included in the delivery.

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Constituent	Measuring range	Repeatability
Alcohol	0.00% ... 15.00%	± 0.02%
Extract	0.00% ... 10.00%	± 0.04%
App. extract	0.00% ... 10.00%	± 0.04%
Original wort	0.00% ... 20.00%	± 0.04%
Density	0.95 bis 1.05 g/cm ³	± 0.0002 g/cm ³
Osm. pressure	Calculated value	—

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Ordering data:

Description: **FermentoFlash** incl. Recording printer

Article-No: 3572

Technical data:

Dimensions (WxHxD): 30 x 24 x 33 cm

Weight: approx. 5 kg

Connected loads: 230 V / 115 V (50/60 Hz) 120 VA